

Morinaga Food Allergen Lateral Flow
Rapid Test Pro II for Gluten

Rev. Jan. 2022

Lot number from #PG001 -

Group	Ingredient	Test Result	Group	Ingredient	Test Result	Group	Ingredient	Test Result
Cereals	Rice powder	-	Sea foods	Salmon	-	Coffee·Cocoa·Tea	Coffee bean	-
	Polished rice	-		Mackerel	-		Cocoa powder	-
	Sprouted brown rice	-		Squid	-		Green tea	-
	Glutinous rice	-		Octopus	-	Seaweeds	Roasted laver	-
	Buckwheat	-		Black tiger prawn	-		Brown seaweed	-
	Japanese barnyard millet	+		Snow crab	-		Kelp 200-fold extract	-
	Foxtail millet	-		Scallop adductor	-		Cumin	+
	Common millet	-		Pollack roe	-	Poppy seeds	-	
	Wheat	/		Salmon roe	-	Black pepper	-	
	Rye	※1		Beef	-	White pepper	-	
	Barley	※1	Chicken	-	Red pepper	-		
	Malt	※1	Pork	-	Wasabi (green horseradish)	-		
	Oat	+	Chicken egg	-	Ginger	-		
	Adlay※2	-	Processed chicken egg	-	Curry powder (spice mix)	-		
	Corn	-	Dried chicken egg	-	Cinnamon	-		
	Corn flour	-	Dairy products	Milk	-	Coriander	-	
	Amaranth	-		Skim milk	-	Turmeric	-	
	Quinoa	-	Green stuff	Tomato	-	Polysaccharide Thickener	Low methoxyl pectin 50-fold extract	-
Sago palm powder	-	Onion		-	Guar gum 1000-fold extract		-	
Yam	-	Spinach		-	Others	Gelatin	-	
Potato starch	-	Fungus	Mushroom	-				
Beans	Soybean		-	Orange	-			
	Red bean (azuki)		-	Kiwi	-			
	Common bean		-	Apple	-			
	Concord pole		-	Peach	-			
	Broad bean		-	Banana	-			
Seeds	Peanut		-	Coconut milk	-			
	Roasted almond	-						
	Roasted cashew nut	-						
	Roasted macadamia nut	-						
	Roasted pistachio nut	-						
	Roasted walnut	-						
	White sesame	-						
Black sesame	-							

Test Result	
-	Negative
+	Positive (reaction positive)

N.T. : Not Tested

※1 These foods have a cross reactivity with this Rapid Test Pro II for Gluten. Please note, if a sample contain a higher concentration of these foods, the test result could be false negative.

※2 The Test Solution may be difficult to flow on a test stick.

These samples were tested by following the instruction of Rapid Test Pro II for Gluten (using Recommended Extraction Method), except Polysaccharide and Thickener samples.

These results can be changed by the difference of kit's lot number or the sample's species or production area.