

Morinaga Food Allergen Lateral Flow
Rapid Test Pro II for Gluten

Rev. Sep. 2022

Lot number from #PG023 -

Group	Ingredient	Test Result	Group	Ingredient	Test Result	Group	Ingredient	Test Result
Cereals	Rice powder	-	Seeds	Peanut	-	Fruits	Orange	-
	Polished rice	-		Roasted almond	-		Kiwi	-
	Sprouted brown rice	-		Roasted cashew nut	-		Apple	-
	Glutinous rice	-		Roasted macadamia nut	-		Peach	-
	Buckwheat	-		Roasted pistachio nut	-		Banana	-
	Japanese barnyard millet	+		Roasted walnut	-		Coconut milk	-
	Foxtail millet	+		Roasted pecan nut	-	Coffee bean	-	
	Common millet	+		White sesame	-	Cocoa powder	-	
	Wheat	+		Black sesame	-	Green tea	-	
	Rye	+		Sea foods	Salmon	-	Roasted laver	-
	Barley	+	Mackerel		-	Hijiki (boiled)	-	
	Malt	+	Squid		-	Brown seaweed (raw)	-	
	Oat	+	Octopus		-	Kelp 200-fold extract	-	
	Adlay	- *	Black tiger prawn		-	Cumin	+	
Corn	-	Snow crab	-		Poppy seeds	-		
Corn flour	-	Scallop adductor	-		Black pepper	-		
Amaranth	-	Pollack roe	-		White pepper	-		
Quinoa	-	Salmon roe	-		Red pepper	-		
Tubers· Starch	Sago palm powder	-	Meats		Beef	-	Ginger	-
	Yam	-		Chicken	-	Cinnamon	-	
	Potato starch	-		Pork	-	Coriander	-	
Beans	Soybean	-	Eggs	Chicken egg	-	Polysaccharide Thickener	Turmeric	-
	Red bean (azuki)	-		Processed chicken egg	-		Low methoxyl pectin 50-fold extract	-
	Common bean	-		Dried chicken egg	-	Guar gum 1000-fold extract	-	
	Concord pole	-	Dairy products	Milk	-	Others	Gelatin	-
	Broad bean	-		Skim milk	-			
Green stuff	Fungus	Tomato	-					
		Onion	-					
		Spinach	-					
Mushroom	-							

Test Result	
-	Negative
+	Positive (reaction positive)

*The Test Solution may be difficult to flow on a test stick.

These samples were tested by following the instruction of Rapid Test Pro II (using Recommended Extraction Method), except Kelp 200-fold extract, Polysaccharide and Thickener samples.

These results can be changed by the difference of kit's lot number or the sample's species or production area.